

MICROLINKS TECHNOLOGY CO.,LTD

# User's Manual

---

## Microscope Installation Manual

MLTC

2014/05/22 Update Ver.2.8A

## Content

<b>1. Please insert CD in CD-Rom .....</b>	<b>3</b>
<b>2. Install and uninstall microscope software.....</b>	<b>3</b>
2.1 Install software.....	4
2.1.1 Setup software .....	4
2.1.2 Software Setup Wizard .....	4
2.1.3 Choose installation route .....	5
2.1.4 Installing.....	5
2.1.5 Install Completed .....	6
2.1.6 Software icon on desktop .....	6
2.2. Uninstall Microscope software .....	7
2.2.1 Choose to uninstall Microscope software.....	7
2.2.2 Confirm to Uninstall .....	7
2.2.3 Uninstalling.....	7
2.2.4 Uninstall Completed.....	7

## 1. Please insert CD in CD-Rom

The package provides setup CD. If the setup CD is missing or damaged, please download setup files from <http://www.vitiny.com>. Please see the following steps to complete the Microscope installation.

- **Step 1: Please insert CD.**
- **Step 2: Install Microscope software. Please refer to section 2.1.**

To uninstall Microscope software in your PC, please follow below 2 steps to uninstall.

- **Step 1: Uninstall Microscope software. Please refer to section 2.2.**

The install window will pop out automatically once CD is inserted.



Fig 1-1 Install window

If the install window does not show, please click Autorun.exe from CD, the same install window will show up, too.



Fig 1-2 Autorun.exe

## 2. Install and uninstall microscope software

Please select the Fig.1-1 to start setup Microscope software program.

## 2.1 Install software

### 2.1.1 Setup software

First, click “Install Microscope software”, next, please select a language, see Fig. 2-1, the initial language will follow up the pc OS system. If you want to cancel setup please click “Cancel” then click “OK”

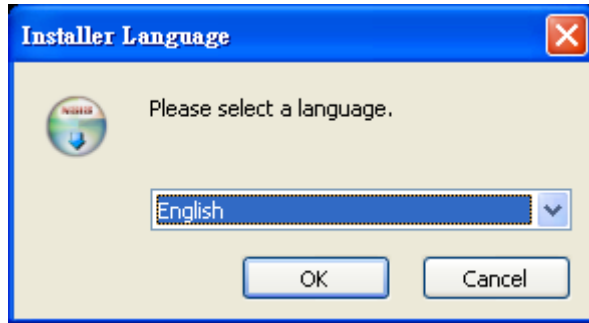


Fig 2-1seup language window

### 2.1.2 Software Setup Wizard

After choosing the language, go to welcome to the Microscope setup wizard window see Fig.2-2, then please click “Next” then continue. If want to cancel please click “Cancel” button. In the following figs, “name” represents Microscope’s software name, “X.XXX.XXXx” represents the software version.

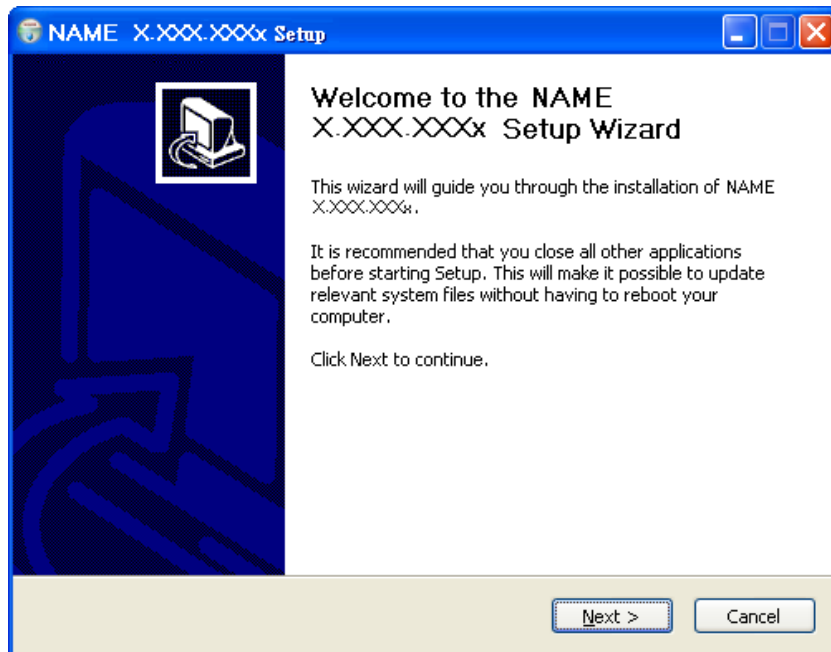


Fig 2-2 welcome to the Microscope setup wizard

### 2.1.3 Choose installation route

Next step please choose the installation route window see Fig.2-3. If want to change file route click “Browse” to change setup path, then please click “install” then start installing.

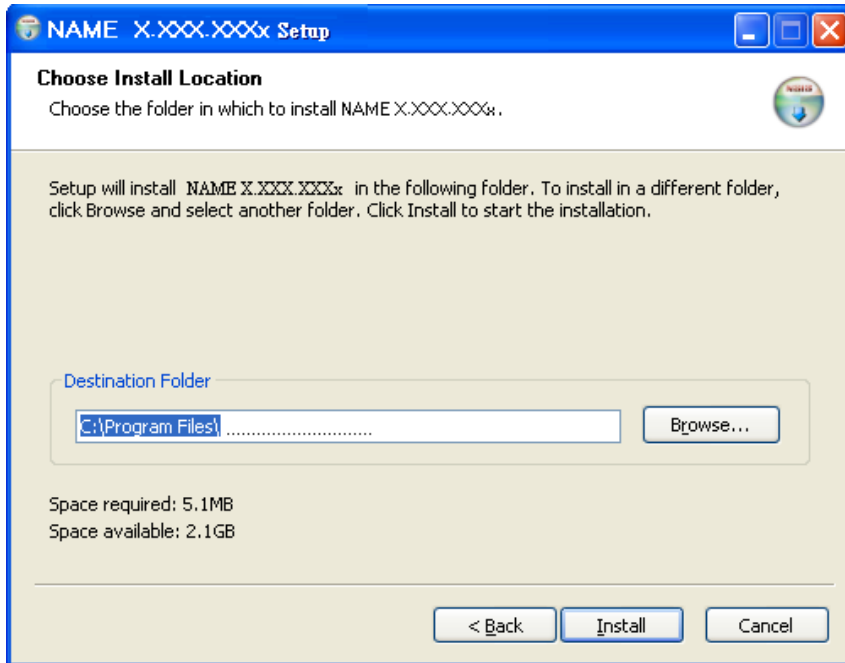


Fig 2-3 choose installation route

### 2.1.4 Installing

Continue install wizard will start installing program ad directory Fig 2-4.

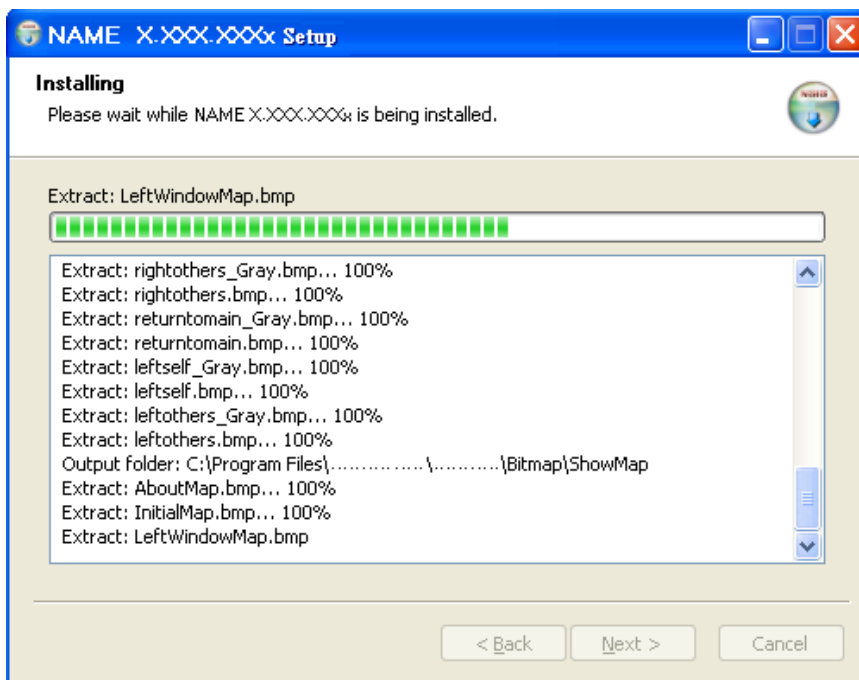


Fig 2-4 installing

### 2.1.5 Install Completed

When the Microscope program complete installation (Fig 2-5), please click “Finish”.

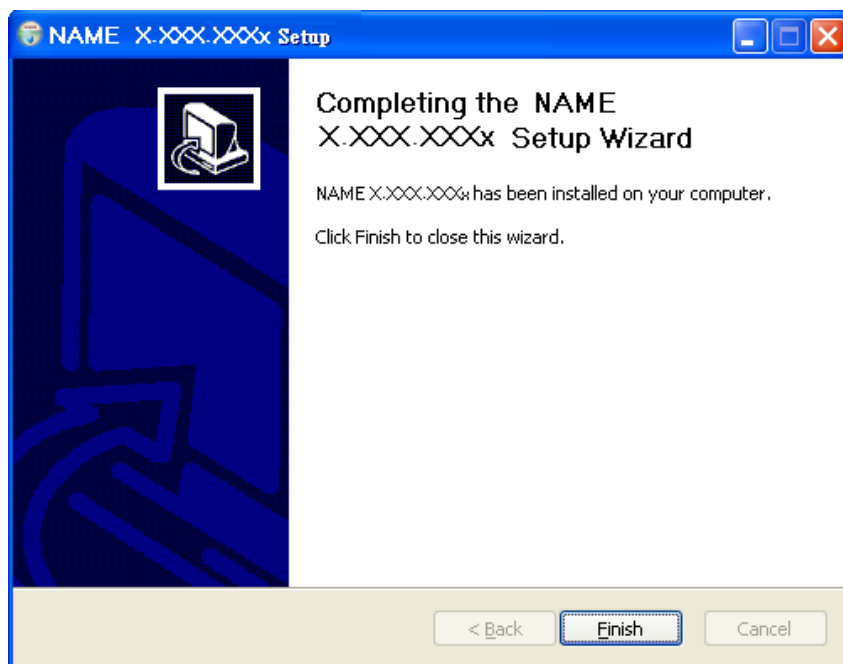


Fig 2-5 Microscope wizard install completed

### 2.1.6 Software icon on desktop

Microscope software shortcut will show on the Desktop as Fig.2-6.



Fig 2-6 Microscope shortcut icon

Microscope shortcut (Fig 2-6) will show on the Desktop. Or go “Start”-> All Programs-> MicroLinks Technology -> Microscope (see Fig 2-7).

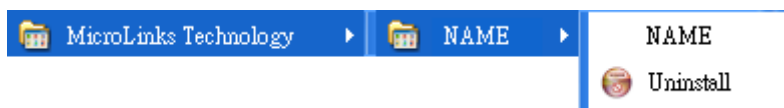


Fig 2-7 Microscope in All Programs

## 2.2. Uninstall Microscope software

### 2.2.1 Choose to uninstall Microscope software

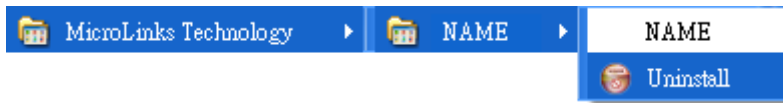


Fig.2-8 Microscope software Uninstall

### 2.2.2 Confirm to Uninstall

Please make sure if you want to completely remove Microscope software. See Fig.2-9

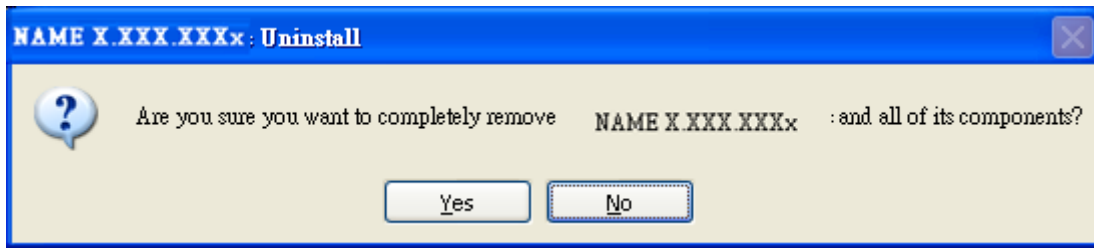


Fig.2-9

### 2.2.3 Uninstalling

In processing of uninstall the Microscope software, see Fig.2-10

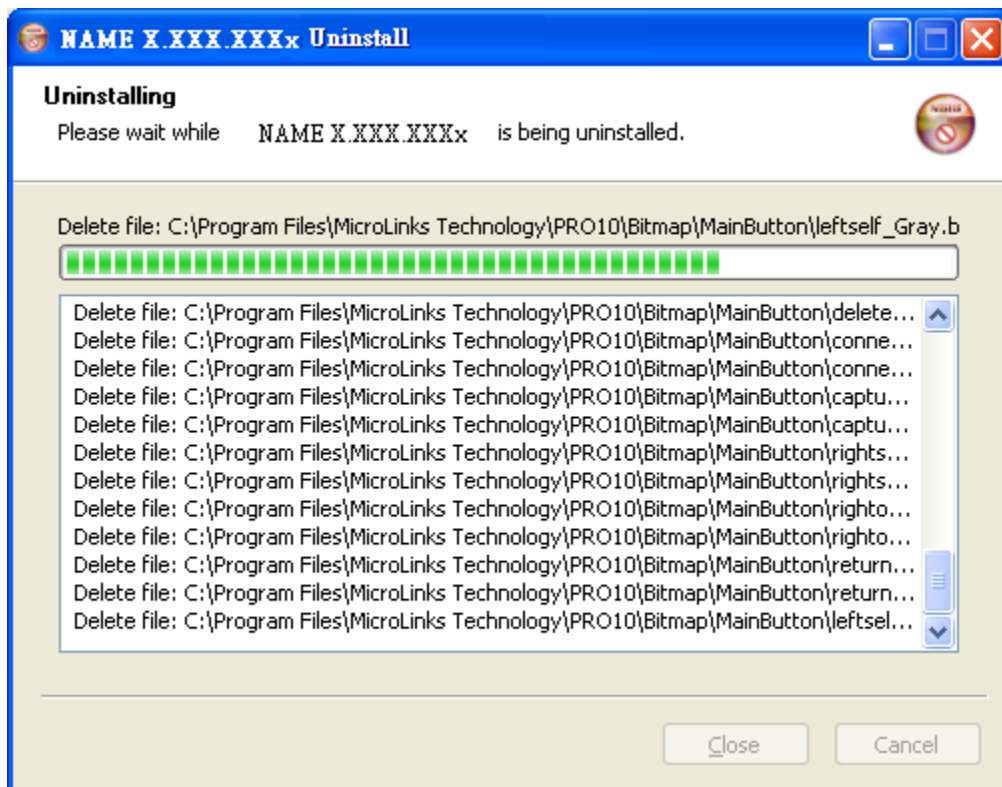


Fig.2-10 In process

### 2.2.4 Uninstall Completed

Microscope software was successfully removed from your PC, please click OK to finish.



Fig. 2-11 Uninstall Completed